


## MC-Media Pad® Coliform test


### Interpretation Guide

The MC-Media Pad® is a convenient method for indicator organism testing. Compliant with international food and beverages standards (AOAC-PTM, MicroVal, ISO 16140), MC-Media Pad® tests streamline your microbial quality control workflow.


#### Accurate results in 4 steps

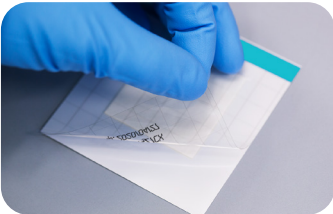


1. Open the cover film diagonally

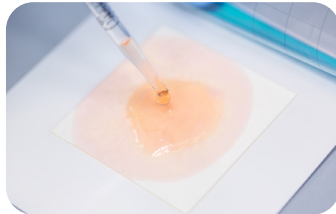

2. Inoculate your sample, it will diffuse automatically

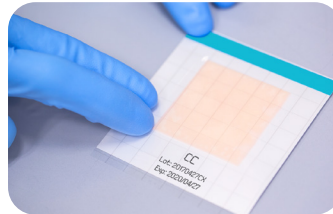

3. Close the cover film and incubate

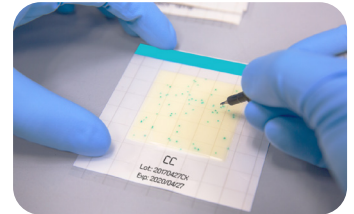

4. Count your results



*Enterobacter aerogenes*  
ATCC® 13048™ 300 CFUs



*Enterobacter aerogenes*  
ATCC® 13048™ > 500 CFU's



*Enterobacter aerogenes*  
ATCC® 13048™ 150 CFUs



*Enterobacter aerogenes*  
ATCC® 13048™ 50 CFUs

With MC-Media Pad®, it is easy to see when a dilution is needed.



Dilution for easier read-out



Dilution factor 1:10



Dilution for easier read-out



Dilution factor 1:10



Ordering information	Product name	Packaging
1.32356.0001	MC-Media Pad® Coliform	100 pcs

Please visit: [SigmaAldrich.com/MC-MediaPad](https://SigmaAldrich.com/MC-MediaPad)

